12 RESPONSIBLE
CONSUMPTION
AND PRODUCTION

Vestfrost Solutions is working towards reaching the UN - Global Sustainable Development Goals by 2030.

The Sustainable Development Goals are the blueprint to achieve a better and more sustainable future for all. In order to implement Goal no 12 "Responsible Consumption and Production", this manual has been printed on recycled paper.

ILR Refrigerator Model VLS 204A/304A/354A/404A/504A AC
Internal Voltage Stabilizer

GB Instructions for use
FR Consignes d'utilisation
ES Instrucciones para el uso
RUS Инструкция по эксплуатации
CHI 使用说明

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WARNING

As the appliance contains flammable refrigerant, as stated on nameplate, it is essential to ensure that the refrigerant pipes are not damaged.

The quantity and type of the refrigerant used in your appliance is indicated on the rating plate.

Standard EN378 specifies that the room in which you install your appliance must have a volume of 1m³ per 8 g of hydrocarbon refrigerant used in the appliances. This is to avoid the formation of flammable gas/air mixtures in the room where the appliance is located in the event of a leak in the refrigerant circuit.

WARNING:

Ventilation openings in the appliance or in built-in structures must be kept clear.

WARNING:

Do not use other mechanical devices or means to accelerate the defrosting process or to remove rime other than those recommended by the manufacturer.

WARNING:

Do not damage the refrigerant system.

WARNING:

The appliance may not be exposed to rain.

WARNING:

This appliance is not intended for use by persons (including children) with reduced physical, sensory or mental capabilities, or lack experience and knowledge, unless they have been given supervision or instructions concerning use of the appliance by a person responsible for their safety. Children should be supervised to ensure that they do not play with the appliance.

WARNING:

Children must not play with, on, or around the appliance.

WARNING:

Children must not clean the appliance or carry out general maintenance unless they are at least 8 years old and are being supervised.

WARNING:

Always, keep the keys in a separate place and out of reach of children.

WARNING:

Before servicing or cleaning the appliance, switch off circuit breaker.

**WARNING:**

Danger risk of fire or explosion. Flammable refrigerant used, as stated on nameplate. To be repaired only by trained personnel.

WARNING:

Do not store explosive substances such as aerosol cans with a flammable propellant in this appliance.

WARNING:

When positioning the appliance, ensure the power cord is not trapped or damaged.

WARNING:

Do not locate multiple portable socket-outlets or portable power supplies at the rear of the appliance.

WARNING:

Appliance use flammable insulation blowing gas. For information about safe disposal, please contact your local disposal service. See section for Disposal.

Contents

WARNING2

Get to know your Refrigerator5

Electrical connection.....6

Voltage stabilizers.....6

HIGHLY IMPORTANT!.....8

Placing the Fridge-tag ®.....9

General information10

Operation and function11

Loading the appliance12

Maintenance and cleaning.....13

Condensation water drainage.....14


Trouble shooting.....15

Warranty, spare parts and service16

Disposal.....17

PQS Code	Model	PQS Performance specifications Specification reference:	PQS Independent type-testing protocol Product verification protocol:
E003/109	VLS 204A AC	E003/RF03.4	E003/RF03-VP.3
E003/110	VLS 304A AC	E003/RF03.4	E003/RF03-VP.3
E003/111	VLS 354A AC	E003/RF03.4	E003/RF03-VP.3
E003/112	VLS 404A AC	E003/RF03.4	E003/RF03-VP.3
E003/113	VLS 504A AC	E003/RF03.4	E003/RF03-VP.3

Get to know your Refrigerator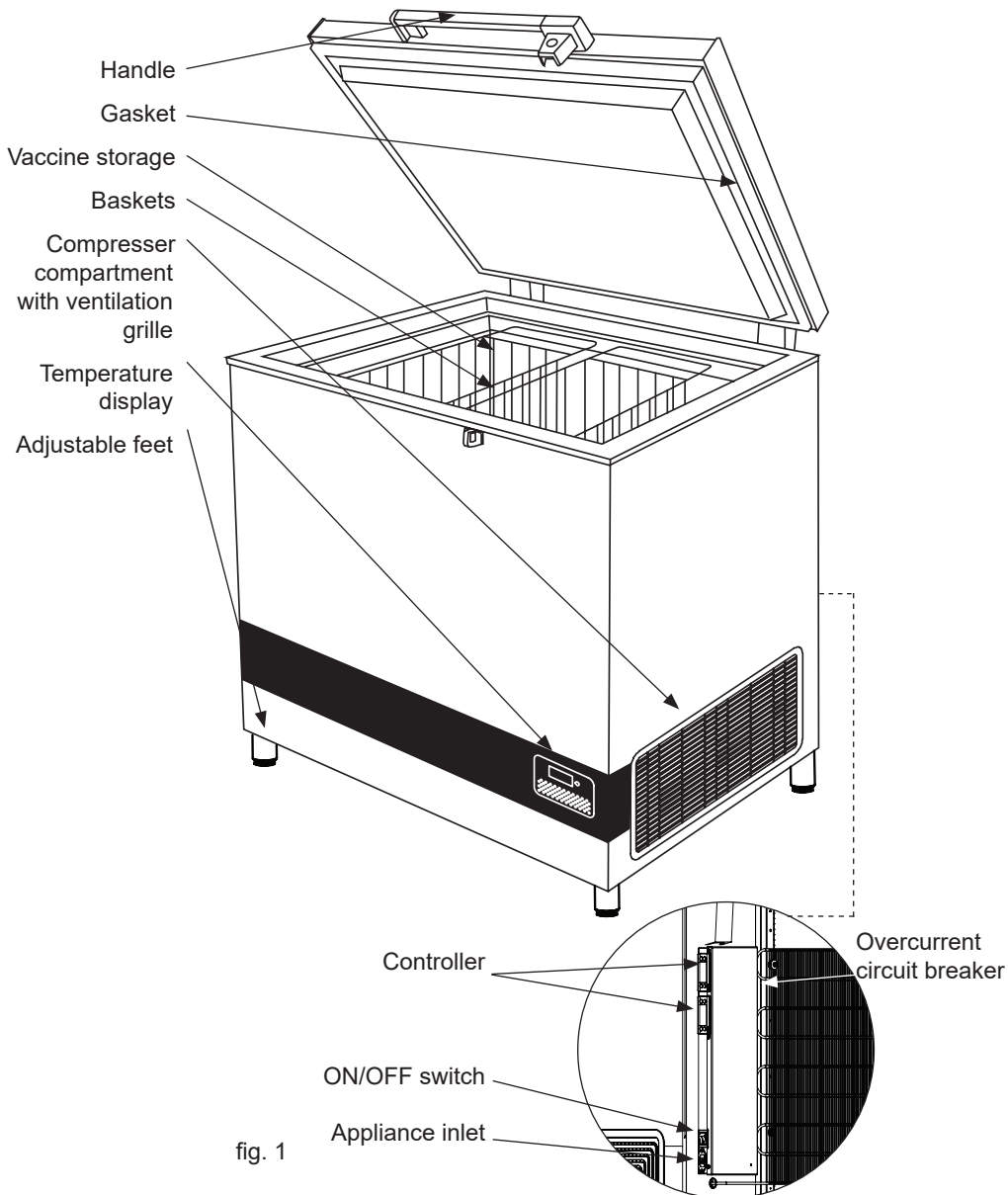

Electrical connection

This device complies with relevant EU directives including Low Voltage Directive 2014/35/EU and Electromagnetic Compatibility Directive 2014/30/EU

The socket should be freely accessible.

Connect the appliance only to 220-240V 50/60Hz mains supply via a correctly installed earthed socket.

The socket must be fused with a 10-13 A fuse.

If the appliance is to be operated in a non-European country, check on the rating plate whether the indicated voltage and current correspond to the values of your mains supply.

Information regarding voltage, current or power are given on the rating plate

Check that the plug fits your type of socket. If it does not fit, procure an appliance cord with the correct plug or have a qualified electrician to fit the plug according to country standards. This appliance must be earthed.

Do not plug several appliances and the refrigerator into the same multiple power strip. The refrigerator should always be plugged into its own individual electrical outlet which has a voltage rating that matches the voltage listed on the refrigerator's rating plate. Ensure that the plug and socket matches and that the appliance is earthed.

Voltage stabilizers

In order to protect the electrical components, the appliance is equipped with an internal voltage stabilizer for extended input range, that will safeguard the equipment against high/low mains voltage, high/low spikes and surges protection in line with WHO/PQS E007/VS01-VP.5

The internal extended stabilizer monitors the mains voltage continuously. If the voltage rises or drops, the stabilizer will stabilize the output to ensure the voltage reaching the appliance remains constant at 230V (+/-2%), within the operating range of the unit. If the input voltage falls below 110V or rises above 285V, the stabilizer will disconnect the output, thereby protecting the load. Once the mains voltage returns within acceptable limits, the stabilizer will reconnect the output following a startup delay.

Delay in restoring supply:

When under or over-voltage cut-out has occurred and the input voltage has returned to the operating range, the output is automatically restored after the factory set delay time delay (3-3:30 minutes random) when the mains voltage again is within the recommended range and stable.

Standard operation instructions how to connect your appliance:

1. Connect the appliance cable to the junction box
2. Ensure that the switch on the appliance is turned OFF
3. Plug the appliance plug into the wall mains socket
4. Turn the power ON at the wall main socket
5. Turn the switch on the appliance ON
6. The internal stabilizer has a built-in AVS (Automatic Voltage Switcher). This will ensure that the load is not connected immediately. The delay will ensure that the mains is good before connecting the load and protects the equipment from rapid switching on and off.
7. Once the waiting period (3-4 min.) has passed, the stabilizer will release power to the appliance, and the cooling system will turn on

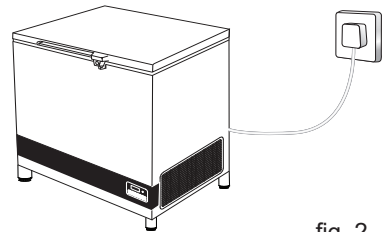
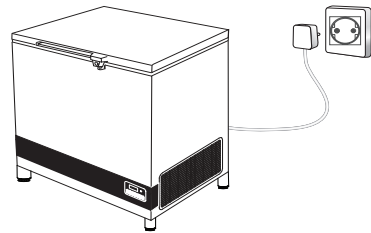
fig. 2

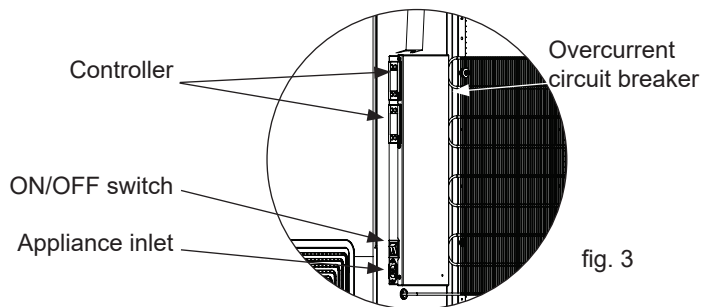
fig. 3

WARNING!

HIGHLY IMPORTANT!

- Make sure NEVER to place any commodities up against the sensor cover.

Eg. Water-pack

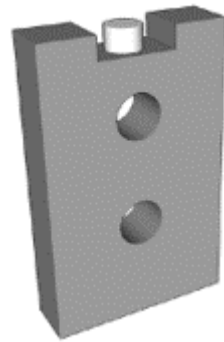
fig. 4

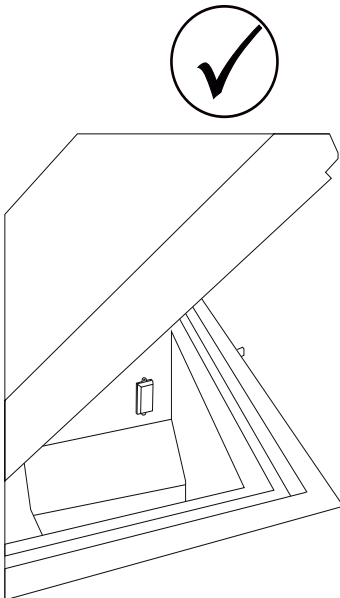
fig. 5

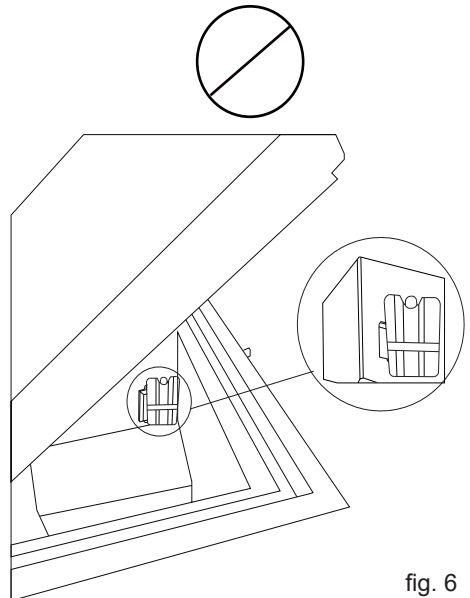
fig. 6

Placing the Fridge-tag[®]

The activated Fridge-tag must be placed immediately in its predetermined location. It is recommended and important to place the device in the centre of the refrigerator for an optimal temperature observation.

External sensor

Two hours before activating the device the external sensor must be placed in its predetermined location. It is recommended and important to place the external sensor in the centre of the refrigerator for an optimal temperature observation and to avoid any incorrect measurements when starting the device

fig. 7

Fridge-tag Support

General information

The VLS A AC is an ice-lined vaccine refrigerator. It has one compartment for vaccine storage (+2 to +8°C) equipped with baskets that are recommended to be used.

Power supply

The VLS A AC line is equipped with a heavy-duty compressor that can operate at voltages 22% below nominal voltage. See the rating plate, which is placed at the rear side of the appliance.

Unpacking

1. Unpack the appliance and check that it has not been damaged. If you observe any damage, inform your supervisor.
2. Check the rating for correct voltage and frequency.
3. Open and remove all internal packing material.

Adjustable feet

Level the appliance by screwing the adjustable feet up or down.

If the appliance is to be placed on a soft surface, e.g. floorboards or a carpet, it is best to recheck whether the appliance is still level after a period of time as the underlying surface may give under the weight of the appliance.

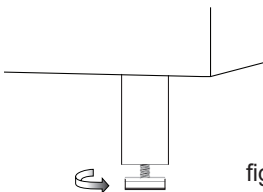
fig. 8

Location

The appliance must be placed in a well ventilated room, not in direct sunlight and away from other heating sources.

When installing the appliance please ensure that the floor is level. Please see figure 9 for further directions.

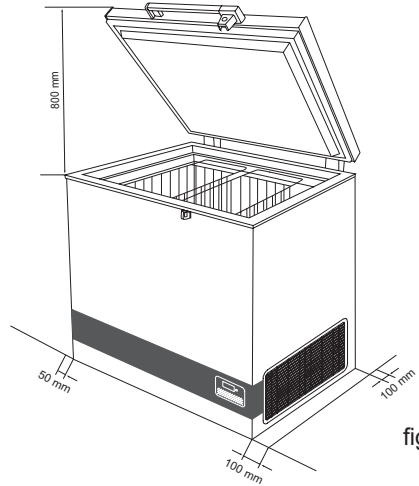
fig. 9

There must be at least 30 mm clearance between the base of the appliance and the floor.

IMPORTANT! There must be free access to the ventilation grille (fig.10).

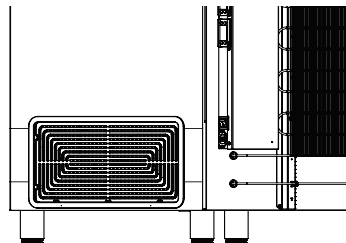
fig. 10

Operation and function

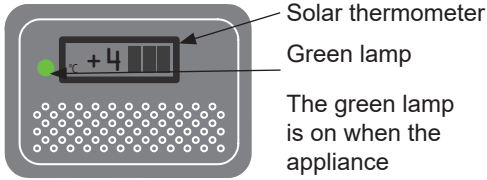
fig. 11 The green lamp is on when the appliance is connected to the power supply.

Starting procedure

Connect the mains lead to the power supply. The green lamp must light (fig. 11) to indicate that the appliance is operating. The electronic thermostat makes a self test, 20 seconds before the compressor starts.

Temperature control

It is recommended to check the inside temperature with an accurate thermometer twice a day. The inside temperature must be checked regularly according to WHO's standards and specifications.

Cool down of the appliance

Before the appliance is loaded with vaccines the ice-lining must be frozen. To ensure that the ice-lining is frozen do the following:

1. Place a thermometer in the top basket.
2. Let the refrigerator run for at least 48 Hours before loading vaccine.
3. Check the temperature in the vaccine compartment (must be between +2° and +8°C).

The temperature in the vaccine compartment must always be monitored on the thermometer and be within the range +2° to +8°C. Due to tolerance of the thermostat you should always control the temperature during cool down.

Controller

The appliance has a controller that controls the temperature in the vaccine compartment. The controller is factory adjusted to a default setpoint, what in most ambient environments will keep the vaccine compartment within desired thresholds of 2°-8°C.

If the cooler is not used for a longer period of time, it is recommended to use the

ON/OFF switch  to turn the cooler OFF. This disconnects the power from mains

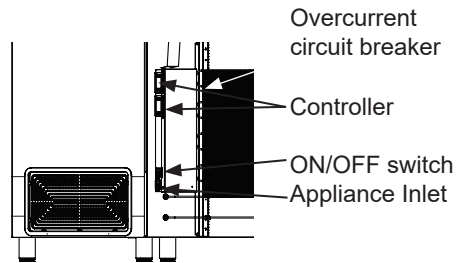
fig. 12

In case cooler is turned OFF, make sure the vaccine compartment is cleaned and dried thoroughly afterward's, this to prevent problems with odors or mold. To prevent odors taking hold it is highly recommended to keep the lid slightly open.

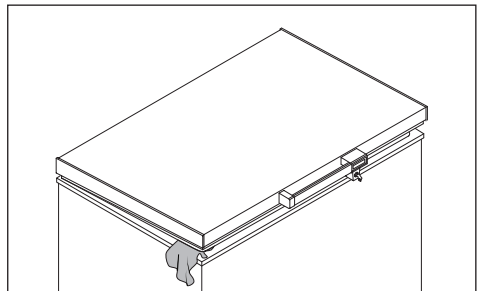
fig. 13

Loading the appliance

Loading vaccines

When the temperature in the vaccine compartment has stabilized, i.e. and the temperature is between $+2^{\circ}$ and $+8^{\circ}\text{C}$ and the compressor stops and starts, vaccines can be loaded. The vaccines should be placed and arranged in the baskets.

To ensure air-circulation and prevent too low vaccine temperature, direct contact to the inside walls must be avoided. The Vaccine load in the baskets shall not be above top of the baskets.(fig. 15)

Vaccine storage

Below points are recommended vaccine storage advice outlined by Unicef

- Keep the vaccine boxes containing the vaccines in neat rows.
- Different vaccines should be kept separately to facilitate easy identification.
- Keep about, 2 cm. space between boxes of vaccines for circulation of air.
- Store Freeze sensitive vaccines (DPT, TT, IPV, Penta and Hep. B) away from the bottom of the ILR to avoid freezing. Always keep the vaccines in the basket provided in the ILR. OPV, RVV BCG, JE and Measles vaccine to be stored at bottom of basket of the ILR.
- Vaccine should be stored as per their heat and cold sensitivity.

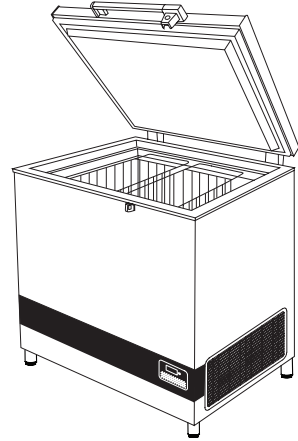
fig. 14

Do not load the vaccine above top of the baskets.

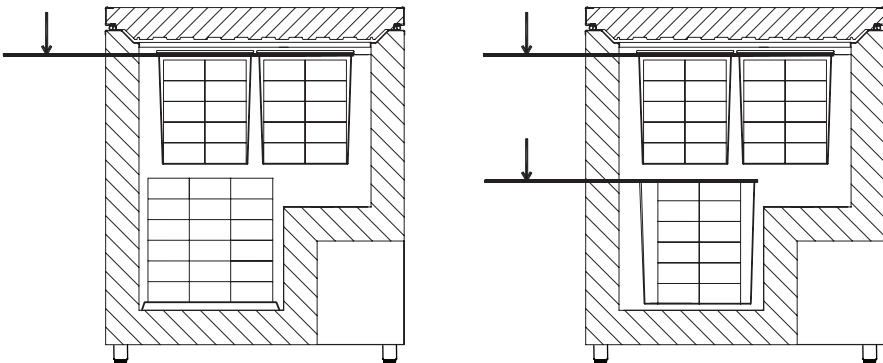
fig. 15

Maintenance and cleaning

Daily maintenance:

The temperature in the vaccine compartment must always be monitored on the thermometer every day. Please note that the vaccine in the appliance must not freeze.

Weekly maintenance:

During normal use, water can accumulate at the bottom of the appliance; remove with a cloth or through the drainage hole. Wipe off water droplets on the inside wall at the same time. Check if lid gasket is sealing tight to the top frame when the lid is closed. A tight sealing lid reduces accumulation of water and formation of rime significantly.

Monthly maintenance:

Clean the grille on the right side of the refrigerator once each month. Clean vaccine compartment with lukewarm water and mild detergent. Clean outside of refrigerator with lukewarm water and mild detergent.

Yearly maintenance:

Electric connections and components are to be checked and cleaned once a year or more if necessary.

Cleaning

Disconnect the power supply before cleaning. The best way to clean the appliance is by using luke warm water with a small amount of unscented detergent. Never use cleaning agents that scour. Use a soft cloth. Rinse with clean water and dry thoroughly. It is important to prevent water from running into the control panel.

The gasket around the lid must be cleaned regularly to prevent discolouration and prolong service life. Use clean water. After cleaning the sealing strip, check that it continues to provide a tight seal.

If the appliance is not being used for any period of time, switch off the appliance, disconnect the power supply, empty the appliance, clean the inside, and leave the lid open to allow air circulation and prevent bad odor.

Condensation water drainage

Due to humidity in the air combined with cold surfaces inside the vaccine compartment, it should be expected that condensation will form on the sides of the inner lining. The condensation will be collected at the bottom, from where it regularly needs to be drained. If the bottom of the vaccine compartment is covered with condensation water, drainage/drying should be initiated.

Condensation increases if:

- Equipment is opened too frequently.
- Lid not closing properly.
- Lid gasket is defective.
- High level of humidity.

Drainage of condensation water is recommended to be performed in the morning between 7am to 9am when the ambient temperatures are expected to be lowest. Proceed as follows:

IMPORTANT!

Before condensation drainage vaccines must be moved to another working ILR or cold box with conditioned icepacks.

1. Disconnect the power supply by switching off the ON/OFF power button or disconnect plug from power socket.
2. Open the drainage plug, placed inside the vaccine compartment.
3. Place a tray under the condensation water drain.

4. When all the condensation and water droplets on the lining has been removed carefully, clean the inside cabinet, as well as lid, lid gasket, and outside of the cabinet, using water with a mild, perfume-free detergent. Dry the vaccine compartment with a soft cloth, and make sure, no moisture is remaining after cleaning.
5. Reinsert the internal drainage plug, and check lid gasket is not faulty.
6. Reconnect the power supply by the ON/OFF power button and/or reconnect power plug.
7. When appliance has regained required safe temperatures for storing vaccines 2° - 8 °C, place the vaccines in neat rows with space.

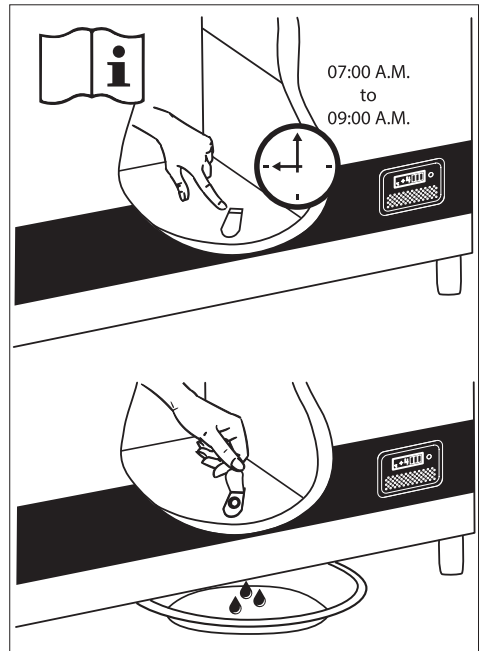
fig. 16

Trouble shooting

Fault	Possible cause	Remedy
Compressor is not running	Be patient, it is most likely that the compressor will start within a few minutes.	If this is not the case, check the following: - Check that power is connected - Check the fuse and replace it if necessary. - If the above is OK, call technical supervisor.
Compressor is running, and the temperature is too high	The ventilation grille is blocked. The lid is not closed properly. The temperature in the room in which the appliance is installed is too high.	Ensure unhindered air circulation. Ensure that the lid is closed properly. Shield the appliance against direct sun light and ensure more ventilation to the room.
No temperature is displayed	The micro processor control unit is broken Power not connected	Change the micro processor control unit. Check power outlet - Check power plug is connected.

Warranty, spare parts and service

Warranty disclaimer

Faults and damage caused directly or indirectly by incorrect operation, misuse, insufficient maintenance, incorrect building, installation or mains connections. Fire, accident, lightning, voltage variation or other electrical interference, including defective fuses or faults in mains installations are not covered by the warranty.

Repairs performed by others than approved service centres and any other faults and damage that the manufacturer can substantiate are caused by reasons other than manufacturing or material faults are not covered by the warranty.

Please note that changes to the construction of the appliance or changes to the component equipment of the appliance will invalidate warranty and product liability, and the appliance cannot be used lawfully. The approval stated on rating plate will also be invalidated.

Before calling for technical assistance, please check whether you are able to rectify the fault yourself. If your request for assistance is unwarranted, e.g. if the appliance has failed as a result of a blown fuse or incorrect operation, you will be charged the costs incurred by your call for technical assistance.

Spare parts

When ordering spare parts, please state the type, serial and product numbers of your appliance. This information is given on the rating plate (fig. 15). The rating plate contains various technical information, including type and serial numbers.

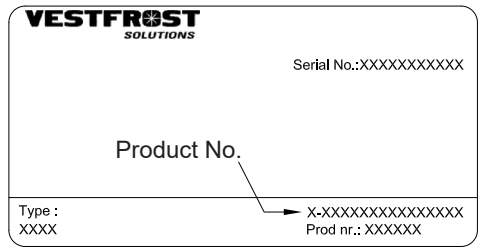
fig. 17

Disposal

Information for Users on Collection and Disposal of Old Equipment and used Batteries

These symbols on the products, packaging, and/or accompanying documents mean that used electrical and electronic products and batteries should not be mixed with general household waste. For proper treatment, recovery and recycling of old products and used batteries, please take them to applicable collection points, in accordance with your national legislation and the Directives 2012/19/EU and 2006/66/EC.

By disposing of these products and batteries correctly, you will help to save valuable resources and prevent any potential negative effects on human health and the environment which could otherwise arise from inappropriate waste handling.

For more information about collection and recycling of old products and batteries, please contact your local municipality, your waste disposal service or the point of sale where you purchased the items.

Penalties may be applicable for incorrect disposal of this waste, in accordance with national legislation.

For business users in the European Union.

If you wish to discard electrical and electronic equipment, please contact your dealer or supplier for further information.

Information on Disposal in other Countries outside the European Union

These symbols are only valid in the European Union. If you wish to discard this product, please contact your local authorities or dealer and ask for the correct method of disposal.

Note for the battery symbol:

This symbol might be used in combination with a chemical symbol. In this case it complies with the requirement set by the Directive for the chemical involved.

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