

Check 3 - Functionality test procedure - Vestfrost CCE AC appliances.

Check	AC	Outcome
1	Wait for the delay set off voltage stabilizer 3-6 min	Yes No
2	Green power light is ON	Yes No
3	Check that controller display is ON	
4	Feel by hand if the compressor is running after 5 minutes of power ON	Yes No
5	Feel by hand if the compressor discharge tube is heating up	Yes No
6	Activate and Note temperature on data logger (Picture to be captured in Vestfrost APP)	°C
7	Feel by hand if the inner lining/evaporator plate is cooling after 1 hour	Yes No
8	After 2 hours Note temperature on data logger (Picture to be captured in Vestfrost APP)	°C

Check 5 - Functionality test procedure - Vestfrost CCE SDD appliances.

Check	SDD	Outcome
1	Green power light is ON	Yes No
2	Check that controller display is ON	Yes No
3	Feel by hand if the compressor is running after 5 minutes of power ON	Yes No
4	Feel by hand if the compressor discharge tube is heating up	Yes No
5	Activate and Note temperature on data logger (Picture to be captured in Vestfrost APP)	°C
6	Feel by hand if the inner lining/evaporator plate is cooling after 1 hour	Yes No
7	After 2 hours Note temperature on data logger (Picture to be captured in Vestfrost APP)	